


- **Highly versatile voltage stimulator for use with all DMT Myograph Systems**
- **Optimized for striated and smooth muscle stimulation**
- **Stimulator is controlled by the included MyoPULSE software delivered with the stimulator's**
- **Easy to connect to a computer**
- **Easy to program and run single pulse, continuous frequency and complicated train stimulation protocols**
- **No need for additional amplifiers**
- **Built-in input and output trigger functions**

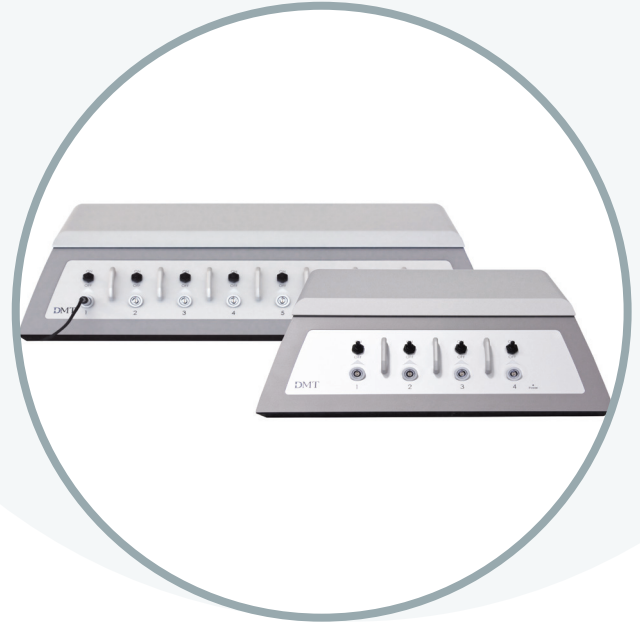

The CS4 and CS8 are 4- and 8-channel Stimulators, respectively. The CS4/CS8 stimulators combine a user-friendly interface with advanced electrical stimulation features required in electro physiological experiments. The CS4 and CS8 are modular, highly versatile voltage stimulators suitable for use with all DMT Myograph Systems.

The CS4 and CS8 have been optimized for striated and smooth muscle stimulation and are ideal for producing controlled perivascular or tissue-embedded neurotransmitter release. The CS4/CS8 Stimulators provide a variety of stimulation modes such as Single Pulse, Continuous Frequency and Train Pulse. The latter is particularly useful when doing fatigue studies in striated muscle and evaluating vascular-neuromuscular junction in smooth muscle

The CS4 and CS8 stimulator are controlled by the MyoPULSE software installed on a PC. MyoPULSE is a flexible software capable of programming simple voltage single pulses and very complicated voltage trains stimulation protocols. All protocol parameters are easily entered and saved in MyoPULSE and can be applied to one or any combination of up to four or eight channel on the CS4 or CS8 stimulator, respectively. All protocols made in the MyoPULSE software can easily be accessed and retrieved.

The CS4 and CS8 stimulators have built-in Input and Output Trigger functions, physically located on the back of the stimulators and controlled by the MyoPULSE software. The Input and Output Trigger functions can be used to trigger external connected equipment e.g. the motor function on a 840MD MyoDYNAMICS Myograph or input trigger to start a stimulation protocol on the CS4 or CS8.



STIMULATOR - CS4/CS8

SPECIFICATIONS:

Output voltage	0-31 Volt
Output	Mono/Bipolar
Stimulus width	0.1-25 m sec.
Max. frequency	2500Hz
Number of pulse in one train	1-32000
Number of trains in one group	1-32000
Number of group	1-32000
Trigger pulse delay	0-3600 sec.
Voltage supply	100-240 Volt 50/60Hz to 48 Volt DC 70 Watt power adapter

